



Trace element solution SL-4 (from medium 14)

Na ₂ -EDTA	0.50	g
FeSO ₄ x 7 H ₂ O	0.20	g
ZnSO ₄ x 7 H ₂ O	0.10	g
MnCl ₂ x 4 H ₂ O	0.03	g
H ₃ BO ₃	0.30	g
CoCl ₂ x 6 H ₂ O	0.20	g
CuCl ₂ x 2 H ₂ O	0.01	g
NiCl ₂ x 6 H ₂ O	0.02	g
Na ₂ MoO ₄ x 2 H ₂ O	0.03	g
Distilled water	1000.00	ml

First dissolve EDTA in distilled water and adjust pH to 7.0 using 2 N NaOH; then add other compounds.