

## 65: GYM STREPTOMYCES MEDIUM

This recipe contains strain-specific modifications for *Ktedonobacter racemifer* DSM 44963 \*

Final pH: \* 5.5

Final volume: 1000 ml

Glucose	4.00	g
Yeast extract	4.00	g
Malt extract	10.00	g
CaCO <sub>3</sub> , if required	2.00	g
Agar	20.00	g
Distilled water	1000.00	ml

Adjust pH to 7.2 before adding agar. Delete CaCO<sub>3</sub> if liquid medium is used.

\* Adjust pH of final medium to 5.5.